



Bacdown[®] Antimicrobial Skin Wipe Technical Data Sheet

- BRIEF DESCRIPTION -

Bacdown Antimicrobial Skin Wipes are for professional laboratory & healthcare use where soap & water are not available. Bacdown Antimicrobial Skin Wipes combine soothing emollients with a broad spectrum antimicrobial agent to provide a gentle wipe that cleans and disinfects skin without drying. Bacdown Antimicrobial Skin Wipes are designed to help minimize infection and contamination risk in laboratory, healthcare, production line, physician, dentist, and veterinarian work places. Bacdown Antimicrobial Skin Wipes meet OSHA criteria for antiseptic hand washing.

- USE PROTOCOL -

Open packet and unfold skin wipe. Wipe hands and fingers for 15-20 seconds to thoroughly wet skin. Repeat whenever necessary.

- ACTIVE INGREDIENTS -

PCMX (Parachlorometaxylene).....0.5%
SD Alcohol 40.....40.0%

- ANTIMICROBIAL ACTIVITY -

Bacdown Antimicrobial Skin Wipes are effective against gram-positive bacteria, gram negative bacteria and fungi. In tests against the 13 organisms listed below, Bacdown Antimicrobial Skin Wipes showed bactericidal, bacteriostatic, fungicidal, and fungistatic activity.

<i>Staphylococcus aureus</i>	<i>Klebsiella pneumoniae</i>
<i>Streptococcus pyogenes</i>	<i>Serratia marcescens</i>
<i>Streptococcus pneumoniae</i>	<i>Candida albicans</i>
<i>Proteus mirabilis</i>	<i>Escherchia coli</i>
<i>Trichophyton mentagrophytes</i>	<i>Salmonella typhi</i>
<i>Pseudomonas aeruginosa</i>	
<i>Methicillin Resistant Staph. aureus</i> (MRSA)	
<i>Vancomycin Resistant Enterococcus faecalis</i> (VRE)	

- SHELF LIFE AND STORAGE -

Bacdown Antimicrobial Skin Wipes have a two-year shelf life when stored at room temperature. The expiration date is printed on each individual wipe's packaging. The wipes are non-flammable.

- OSHA RULES AND REGULATIONS -

When provision of hand washing facilities is not feasible, the employer shall provide either appropriate antiseptic hand cleanser in conjunction with clean cloth/paper towels or antiseptic towelettes. When antiseptic hand cleansers or towelettes are used, hands shall be washed with soap and running water as soon as feasible (Federal Register /Vol. 58, No. 235/Friday, December 6, 1991/Rules and Regulations).

- PRECAUTIONARY STATEMENT -

For external use only. Keep out of eyes and avoid contact with meninges. If Bacdown Antimicrobial Skin Wipe liquid should get into eyes, rinse eyes promptly and thoroughly with water.

- PACKAGING -

50 wipes per pack / 20 packs per case
Dimensions of each wipe are 5 inches by 8 inches.

- CATALOG NUMBER -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	20 packs/cs	7003

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

©Decon Labs, Inc.

Revised 11/16/06