



CiDecon® Plus Wipes

Surface Disinfectant Wipes

Technical Data Sheet

EPA Reg. No. 8383-7-56753
CANADA DIN # 02203294

- BRIEF DESCRIPTION -

CiDecon Plus Wipes are disinfectant wipes in a convenient pop-up canister. An effective disinfectant and active deodorant for laboratories, hospitals, clinics, medical, dental and veterinary offices, industrial clean rooms, homes, nursing homes, ambulances, etc. For use on hard inanimate surfaces, they provide broad range efficacy against bacteria, virus and fungi. Proven effective against tuberculosis, HIV-1 and many other pathogenic organisms. CiDecon Plus Wipes are safe, effective and meet OSHA bloodborne pathogen standard.

- DIRECTIONS FOR DISPENSING -

Pull up tab and remove cover from container. and pull up the wipe corner from the center. Twist the corner into a point and insert about one inch of the wipe through the X hole on the underside of the cover. Pull the wipe completely through and replace cover onto container. Pull out first wipe at a 90° angle. Remaining wipes feed automatically. Always reseal the hinged cap after each use to ensure moisture retention.

- USE PROTOCOL -

It is a violation of U.S. Federal Law to use this product in a manner inconsistent with its labeling. PRECLEAN surface to remove soil and filth. Wipe dry.

- ACTIVE INGREDIENTS -

Phenol 1.56 %
Sodium Phenate 0.06 %
Inert Ingredients 98.38 %

- EFFICACY -

CiDecon Plus Wipes act effectively in three minutes for hospital disinfection. CiDecon Plus Wipes also provide a continuous residual bacteriostatic activity for six months.

CiDecon Plus Wipes are germicidal when used as directed.

PASSES AOAC EFFICACY STANDARDS FOR INSTITUTIONAL AND HOSPITAL DISINFECTION

- Kills HIV-1 (AIDS virus) on hard, nonporous surfaces in one minute.
- Kills *Salmonella choleraesuis*, *Staphylococcus aureus*, *Pseudomonas aeruginosa* and *Trichophyton mentagrophytes* (Athlete's Foot fungus) on hard nonporous surfaces in 3 minutes.
- Kills vegetative organisms including *Streptococcus pyogenes*, *Streptococcus salivarius*, *Escherichia coli*, *Herpes simplex* types 1/F and 2/G (oral, ocular and genital), *Influenza A₂* (Japan 305/57 Asian strain) *Vaccinia*, *Canine paravirus*, *Cytomegalovirus*, *Coronavirus* and *Polio type 1* viruses on hard nonporous inanimate surfaces and *Mycobacterium tuberculosis* at 20° C/68° F or above in 10 minutes.

- STORAGE AND DISPOSAL -

Store in cool area. Do not reuse wipes. Dispose of used wipe and empty dispenser in trash. Materials that may contain blood and other body fluids should be autoclaved and disposed of according to federal, state and local regulations for infectious waste.

- PRECAUTIONARY STATEMENT -

Hazard to humans and domestic animals.

CAUTION: MAY BE IRRITATING TO EYES.

If in eyes, immediately flush with water for at 15-20 minutes. Remove contact lenses, if present, after the first 5 minutes, then continue rinsing eyes. Get medical attention. If on skin, wash thoroughly with water. Wash contaminated clothing before reuse. **WASH THOROUGHLY AFTER HANDLING WIPES. KEEP OUT OF REACH OF CHILDREN.**

- PACKAGING -

180 wipes per canister / 12 canisters per case

- CATALOG NUMBER -

<u>Manufactured for:</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	180 wipes	8511

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**
