



CiDecon[®] Q

Germicidal Foaming Cleaner

EPA 11694-34-56753

Technical Data Sheet

- BRIEF DESCRIPTION -

CiDecon Q Germicidal Foaming Cleaner is an effective disinfectant / deodorant that can be used on most hard, non-porous surfaces to clean, disinfect and deodorize all in one step. CiDecon Q has a light lemon fresh scent.

- USE PROTOCOL -

CiDecon Q foam clings to vertical surfaces permitting longer contact time. To clean, remove cap, shake can, spray evenly over the surface and wipe clean. For heavily soiled surfaces allow foam to remain for one minute before wiping.

To disinfect, first clean the surface thoroughly as directed above. Then spray a thin coat of foam, making sure it covers the entire surface to be disinfected. Allow to remain for at least ten minutes, and then wipe off with clean damp cloth. See additional information on product label.

- EFFICACY -

CiDecon Q disinfects the following organisms on hard nonporous inanimate environmental surfaces: *Herpes Simplex Type 1 virus, Herpes Simplex Type 2 (genital) virus, Influenza A₂ (Hong Kong), Human Immunodeficiency virus, HIV-1 (Aids virus) and Human Coronavirus*. This product will also kill the following organisms: *Staphylococcus aureus, Pseudomonas aeruginosa, and Salmonella choleraesuis*.

- STORAGE AND DISPOSAL -

STORAGE. Contents under pressure. Store containers indoors and in a cool dry area away from heat and open flames. Replace cap when not in use.

PESTICIDE DISPOSAL. Do not contaminate water, food or feed by storage or disposal. Wastes resulting from the use of this product may be disposed of on site or at an approved waste disposal facility.

CONTAINER DISPOSAL. Do not puncture or incinerate container.

- PRECAUTIONARY STATEMENT -

HAZARDS TO HUMAN AND DOMESTIC ANIMALS.

Causes substantial but temporary eye injury. Harmful if absorbed through the skin. Avoid breathing spray mist. Do not get in eyes, on skin or on clothing. Wear safety glasses or goggles. Wash thoroughly with soap and water after handling and before eating, drinking, chewing gum, or using tobacco. Remove and wash contaminated clothing before reuse.

-FIRST AID-

If in eyes: Hold eye open and rinse slowly and gently with water for 15-20 minutes. Remove contact lenses, if present, after the first 5 minutes, then continue rinsing eye. Call a poison control center or doctor for treatment advice. If inhaled: move person to fresh air. If person is not breathing, call 911 or an ambulance, then give artificial respiration, preferably mouth to mouth, if possible. Call a poison control center or doctor for treatment advice. If swallowed: Call a poison control center or doctor for treatment advice. Do not induce vomiting unless told to do so by a poison control center or doctor. Do not give anything by mouth to an unconscious person. If on the skin or clothing: Take off contaminated clothing. Rinse skin immediately with plenty of water for 15-20 minutes. Call a poison control center or doctor for treatment advice.

- PACKAGING -
18 oz. / 12 per case

- CATALOG NUMBER-

<u>Manufactured for:</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs	18oz.	8618

For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com

[®]Decon Labs, Inc. Revised: 1/2009