



Contrex® CF is a cation-free and phosphate-free liquid detergent concentrate. Contrex® CF is formulated for use in cleaning of equipment and electronics that are sensitive to sodium and metal cations and also phosphates.

APPLICATIONS:

Contrex® CF is totally rinsable eliminating detergent build-up that interferes with analysis and other critical procedures.

To accelerate cleaning - increase concentration and/or temperature. After cleaning, rinse thoroughly. Cleaning applications include:

- **Electronic components**
- **Solder/Flux removal**
- **Environment lab ware**

**Decon Catalog No. 5801
4 x 1 Gallon**

WORKING SOLUTIONS: Contrex® CF may be diluted with distilled, deionized, or clean potable water and can be used as a soak, scrub, ultrasonic cleaner, or a machine washing. Make a fresh 3% to 5% Contrex® CF working solution (4 to 6oz. per gallon) in deionized or distilled water for best results. Solutions of Contrex® CF are biodegradable and readily disposable. Protective gloves and safety glasses are recommended.

RINSABILITY: Contrex® CF is totally rinsable - however - numerous factors affect rinsability. These factors include water quality, temperature, and volume of water, nature of surface, shape, and size of the vessel.

STORAGE: Store concentrate in a sealed container at ambient temperature. Do not freeze.



For details on Decon's Full Line of Contrex Detergents visit

www.deconlabs.com

Rev. 09/2022