



## Technical Data Sheet

---

### - BRIEF DESCRIPTION -

*Contrex TL effervescent detergent tablets are easy to use with no measuring required. Contrex TL is formulated for use in pipet cleaning, parts cleaning or in any cleaning application assisted by effervescent action.*

### - APPLICATIONS -

To wash pipettes in siphon-type washers: If needed pre-soak pipettes immediately after use in a 1 % Contrex AP, Contrex AL, or Contrex EZ solution. Place pipet inside washer and drop one Contrex TL tablet into the pipet washer. Contrex TL tablet will dissolve (in cold or warm water) and effervesce, producing a detergent solution that passes up and down through the bores and around the pipettes. Commence rinsing after the tablet is dissolved

For critical work, use distilled, deionized or purified water for the final rinse. Cleaning applications include pipettes and tubes in siphon washers.

### - RINSABILITY -

Contrex TL is free rinsing and disposable straight to the drain

### - STORAGE -

Store in a clean, dry environment.

### - PRECAUTIONARY STATEMENTS -

**KEEP OUT OF REACH OF CHILDREN**

#### CAUTION

Use only as directed and read label and safety information before using. Personal protective equipment such as safety glasses and gloves is recommended.

### - PACKAGING -

100 tablets / 6 per case

### - CATALOG NUMBERS-

<u>Item</u>	<u>Cat. No.</u>
100 tablets	5900

---

For Additional Technical Information call  
Decon Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)

---

®Decon Labs, Inc.

Revised 11/16/06