



Technical Data Sheet

- BRIEF DESCRIPTION -

Contrex EZ Enzyme Cleaner is a concentrated, powdered enzymatic detergent. It is formulated for use as an enzymatic detergent for use in manual or soak cleaning of labware and equipment contaminated with protein and fats.

- APPLICATIONS -

Contrex EZ is not for use in machines, product will foam. Safe on most surfaces including glass, plastic and most metals. Cleaning applications include:

- Labware
- Instruments
- Endoscopes
- Medical Devices
- Process equipment
- Food and Dairy surfaces
- Ultra filtration and reverse osmosis filters

- WORKING SOLUTIONS -

Contrex EZ concentrate can be diluted with distilled, deionized or clean potable water. Make up a fresh 1 % Contrex EZ working solution (1.25 oz. per gallon) in deionized or distilled water for best results. Clean by scrubbing, soaking or by ultrasonic cleaning.

- RINSABILITY -

Contrex EZ is totally rinsable – however, numerous factors affect rinsability. These factors include water quality, temperature, and volume of water, nature of surface, shape and size of the vessel.

-STORAGE -

Product must be stored in a clean, dry environment.

- PRECAUTIONARY STATEMENTS -

KEEP OUT OF REACH OF CHILDRREN CAUTION

Contents may cause eye irritation. Avoid contact with eyes.

-FIRST AID-

If in eyes: flush eyes with copious amounts of water for 15 minutes. Seek medical attention if eye irritation persists.

- PACKAGING -

4 lb. / 9 per case
25 lb. / each

- CATALOG NUMBERS-

<u>Item</u>	<u>Cat. No.</u>
4 lbs	5404
25 lbs	5425

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

®Decon Labs, Inc.

Revised 4/02/2012