



Germicidal Foaming Cleaner EPA 11694-34-56753 Technical Data Sheet

- BRIEF DESCRIPTION -

CiDecon Q Germicidal Foaming Cleaner is an effective disinfectant / deodorant that can be used on most hard, non-porous surfaces to clean, disinfect and deodorize all in one step. CiDecon Q has a light lemon fresh scent.

- USE PROTOCOL -

To clean: Remove cap, shake can, spray evenly over the surface and wipe clean. For heavily soiled surfaces allow foam to remain for one minute before wiping.

To disinfect: Clean the surface thoroughly as directed above. Then spray a thin coat of foam, making sure it covers the entire surface to be disinfected. Allow to remain for at least 10 minutes, then wipe off with clean damp cloth. See additional information on product label.

- ACTIVE INGREDIENTS -

n-Alkyl (60% C ₁₄ , 30% C ₁₆ , 5% C ₁₂ , 5% C ₁₈) dimethyl benzyl ammonium chlorides.....	0.1%
n-Alkyl (68% C ₁₂ , 32% C ₁₄) dimethyl ethylbenzyl ammonium chlorides	0.1%
Inert Ingredients	99.8%

- EFFICACY -

CiDecon Q disinfects the following organisms on hard,non-porous, inanimate environmental surfaces: *Herpes Simplex Type 1 virus, Herpes Simplex Type 2 (genital) virus, Influenza A₂ (Hong Kong), Human Immunodeficiency virus, HIV-1 (Aids virus), Staphylococcus aureus, Pseudomonas aeruginosa, Salmonella choleraesuis.* and *Human Coronavirus(HCoV).* Kills Pandemic 2009 H1N1 Inflenza A virus (formerly called swine flu).

Kills HIV-1, on pre-cleaned environmental surfaces/objects previously soiled with blood/body fluids in healthcare setting or other settings in which there is an expected likelihood of soiling of inanimate surfaces/objects with blood or body fluids and in which the surfaces/objects likely to be soiled with blood or body fluids can be associated with the potential for transmission of Human Immunodeficiency Virus Type 1 (HIV-1) (associated with AIDS).

SPECIAL INSTRUCTIONS FOR CLEANING AND DECONTAMINATION AGAINST HIV-1 HBV AND HCV ON SURFACES / OBJECTS SOILED WITH BLOOD/BODY FLUIDS:

PERSONAL PROTECTION: When handling items soiled with blood or body fluids use disposable latex gloves, gowns, masks, or eye coverings.

CLEANING PROCEDURES. Blood and other body fluids must be thoroughly cleaned from surfaces and objects before application of CiDecon Q.

CONTACT TIME. Allow at least 10 minutes contact time with the surfaces/objects for disinfection.

DISPOSAL OF INFECTIOUS MATERIAL: Contaminated waste and blood and other body fluids should be autoclaved and disposed according to local regulations for infectious waste disposal. This product is not to be used as a terminal sterilant/high level disinfectant on any surface or instrument that (1) is introduced directly into the human body, either into or in contact with the bloodstream or normally sterile areas of the body, or (2) contacts intact mucous membranes but which does not ordinarily penetrate the blood barrier or otherwise enter normally sterile areas of the body.

- STORAGE AND DISPOSAL -

STORAGE. Contents under pressure. Store in a cool dry area away from heat and open flames. Exposure to temperatures above 120 °F may cause bursting.

CONTAINER DISPOSAL. Do not puncture or incinerate container. If empty: Place in trash or offer for recycling. If partly filled: Call your local solid waste agency for disposal instructions.

**- PRECAUTIONARY STATEMENTS -
KEEP OUT OF REACH OF CHILDREN.**

WARNING

HAZARDS TO HUMAN AND DOMESTIC ANIMALS. Causes substantial but temporary eye injury. Harmful if absorbed through the skin. Avoid breathing spray mist. Do not get in eyes, on skin or on clothing. Wear safety glasses or goggles. Wash thoroughly with soap and water after handling and before eating, drinking, chewing gum, or using tobacco. Remove and wash contaminated clothing before reuse.

FIRST AID: If in eyes: Hold eye open and rinse slowly and gently with water for 15-20 minutes. Remove contact lenses, if present, after the first 5 minutes, then continue rinsing eye. Call a poison control center or doctor for treatment advice. If inhaled: move person to fresh air. If person is not breathing, call 911 or an ambulance. Call a poison control center or doctor for treatment advice. If swallowed: Call a poison control center or doctor for treatment advice. Do not induce vomiting unless told to do so by a poison control center or doctor.

If on skin or clothing: take off contaminated clothing. Rinse skin immediately with plenty of water for 15-20 minutes.

- PACKAGING -

18 oz. / 12 per case

- CATALOG NUMBER-

<u>Manufactured for:</u>	<u>Item</u>	<u>Cat.</u>
<u>No.</u> Decon Labs	18oz.	8618

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

page 2 of 2

®Decon Labs, Inc. Revised: 12/12/2012