



Surface Disinfectant Wipes

Technical Data Sheet

EPA Reg. No. 55195-4-56753

- BRIEF DESCRIPTION -

CiDecon Wipes are disinfectant wipes in a convenient pop-up canister. For use on hard, non-porous, inanimate surfaces, they provide broad range efficacy against bacteria, virus and fungi. Proven effective against HIV-1 and many other pathogenic organisms. CiDecon Wipes are safe, effective and meet OSHA bloodborne pathogens standard.

- DIRECTIONS FOR DISPENSING -

Remove large snap cover from container and pull up the wipe corner from the center. Twist the corner into a point and thread through the X hole on the underside of the cover. Pull the wipe completely through and replace cover onto container. Pull out first wipe at a 90° angle. Remaining wipes feed automatically. Always reseal the hinged cap after each use to ensure moisture retention.

- USE PROTOCOL -

Thoroughly wet pre-cleaned surface with wipe and allow to air dry. Blood and other bodily fluids must be thoroughly cleaned from surface and objects before application of wipes.

- ACTIVE INGREDIENTS -

Glutaraldehyde.....0.27500%
 o-Phenylphenol.....0.01375%
 p-Tertiary Amylphenol.....0.00275%

- EFFICACY -

When used as directed, CiDecon Wipes kill:
(10 minutes): *Salmonella choleraesuis, Staphylococcus aureus, Methicillin- Resistant Staphylococcus aureus (MRSA), Pseudomonas aeruginosa, Trichophyton mentagrophyte, Human Coronavirus (HCoV).*
(2 minutes): *Coxsackie virus B5a, Adenovirus-2, Poliovirus 1 & 2, Rotavirus SA-11, Rhinovirus, Vaccinia virus, Cytomegalovirus, Respiratory Syncytial Virus.*
(30 seconds) HIV-1 (AIDS virus), Herpes simplex virus 1 & 2, Influenza A2HK
 Kills Pandemic 2009 H1N1 Influenza A virus (formerly called swine flu)

- STORAGE AND DISPOSAL -

Store in cool area. Do not reuse wipes. Dispose of used wipe in trash after use. Do not flush. Materials that may contain blood and other body fluids should be autoclaved and disposed of according to federal, state and local regulations for infectious waste. Non-refillable container. Do not reuse container. Offer empty container for recycling if available or discard in trash.

- PRECAUTIONARY STATEMENT -

Hazards to humans and domestic animals.
CAUTION: MAY BE IRRITATING TO EYES.
 Do not get in eyes, on skin or on clothing. Wear rubber gloves and safety glasses when handling. Repeated dermal contact may result in dermal sensitization in some individuals. Harmful if swallowed. Avoid contamination of food. Avoid reuse of unwashed disinfectant-soaked glove. Avoid prolonged inhalation.

STATEMENT OF PRACTICAL TREATMENT
 If in eyes, immediately flush with water for at least 15 minutes. Get medical attention. If on skin, wash thoroughly with water. Wash contaminated clothing before reuse. **WASH THOROUGHLY AFTER HANDLING WIPES. KEEP OUT OF REACH OF CHILDREN.**

- PACKAGING -

65 wipes per canister / 6 canisters per case

- CATALOG NUMBER -

Manufactured for:	Item	Cat. No.
Decon Labs, Inc.	65 wipes	8510

**For Additional Technical Information call
 Decon Customer Service
 (800) 332-6647, or
 See our website at www.deconlabs.com**

®Decon Labs, Inc. Revised 8/18/2011