



CiDehol[®] 70

Filtered 70% IPA Solution

Technical Data Sheet

- BRIEF DESCRIPTION -

CiDehol 70 is a 70% (v/v) Isopropyl alcohol solution filtered to 0.2 microns. CiDehol 70 can be used to clean surfaces in microbiology labs and tissue culture labs. It is also effective in cleaning surfaces and process equipment in pharmaceutical, biotechnology, and medical device, cosmetic and food production areas. IPA solution evaporates away completely so no rinse is required.

- USE PROTOCOL -

Use in well-ventilated area. Spray or wipe solution on surface to be cleaned until it is completely covered. Allow to air dry. When using trigger spray, point away from eyes and skin. Hold bottle six to eight inches from surface to be treated.

- INGREDIENTS -

70% Isopropyl Alcohol, USP
30% Purified Water, USP

-SHELF LIFE AND STORAGE -

The shelf life for CiDehol 70 is 2 years from the date of manufacture. The expiration date can be found on the product. Do not store at temperatures above 120°F.

- PRECAUTIONARY STATEMENT -

Do not come in contact with flame or extreme heat. Avoid eye and skin contact and ingestion.

- PACKAGING -

16 oz. Trigger Spray / 12 per case
16 oz. Trigger Spray / 12 per case (Color Label)
32 oz. /12 per case
1 Gallon / 4 per case
5 Gallons
55 Gallons

- CATALOG NUMBERS -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	16 oz.	8416
	32 oz.	8432
	1 Gal.	8401
	5 Gal.	8406
	55 Gal.	8405
Color Label	16 oz.	8417

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

®Decon Labs, Inc. Revised 11/11/2011