



# ***Dri-Clean*** **Powdered Detergent**

## **Technical Data Sheet**

---

### **- BRIEF DESCRIPTION -**

*Dri-Clean is a powdered detergent for use in cleaning all labware and equipment. Use Dri-Clean for soak, manual, and ultrasonic cleaning applications. Removes protein, dried blood, and most organic and inorganic contaminants from glassware, plasticware, ceramics, and metal surfaces. Completely free-rinsing and residue-free. Dri-Clean is packaged in a wide-mouth plastic container with a convenient handle. This makes it easier to handle and prevents the absorption of water and caking associated with powdered detergents packaged in milk cartons.*

### **- USE PROTOCOL -**

Dissolve 1/2 oz. (1 Tbsp.) in one gallon of water. For tenacious contaminants, use 1 oz. (2 Tbsp.) or more and dissolve in hot water. Clean by manually scrubbing, soaking, or use ultrasonic bath. For critical applications, clean and rinse with distilled or deionized water.

### **- STORAGE -**

Store in original container with lid tightly sealed. Contamination with water or high humidity may cause caking.

### **- PRECAUTIONARY STATEMENT -**

Contains sodium carbonate. May cause eye and skin irritation. Avoid eye contact or prolonged skin contact. Avoid breathing dust.

**WEAR PROTECTIVE EYEWARE AND GLOVES. KEEP OUT OF REACH OF CHILDREN**

### **FIRST AID:**

Eye Contact: Flush eyes with water for 15 minutes. Get medical attention if irritation persists.

Skin Contact: Rinse with water, remove contaminated clothing.

Ingestion: Give water or milk. Never give anything by mouth to an unconscious person. Get medical attention immediately.

### **- PACKAGING -**

2 kg (4.4 lbs) / 6 per case  
50 kg (110 lbs)

### **- CATALOG NUMBERS -**

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	2 kg (4.4 lbs)	2502
	50 kg (110 lbs)	2501

---

**For Additional Technical Information call  
Decon Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)**

---

®Decon Labs, Inc.

Revised 11/16/06