

# DNase Displace<sup>™</sup>

## Technical Data Sheet

---

### - BRIEF DESCRIPTION -

*DNase Displace removes DNase, and unwanted DNA. It works instantly so there is no waiting. DNase Displace eliminates the time-consuming method of autoclaving. DNase Displace can be used on all surfaces including glass, plastic, metal, and ceramic. DNase Displace is phosphate-free, safely disposable, non-toxic, and totally rinsable for the decontaminated surface.*

### - USE PROCEDURE -

**Use DNase Displace full strength.**

**Method 1:** Thoroughly wet item to be decontaminated with DNase Displace - ensure all surfaces are coated. Swirl for a few seconds, drain, and immediately rinse twice with the DNase-free water.

**Method 2:** Soak item in DNase Displace until needed. Drain and rinse as above.

**Method 3:** Soak an DNase-free wipe in DNase Displace, wipe over surface to be decontaminated. Rinse as above.

**For total radioactive decontamination,** soak item in DNase Displace for 2-12 hours. Rinse as above. Wipe over benchtops and sinks to remove residual contamination.

### - CAUTION -

**Alkaline!** Severe eye irritant, mild skin irritant. Wear eye and skin protection when handling DNase Displace. In case of contact with eyes, immediately flush with water for fifteen minutes and contact a physician immediately.

### - PACKAGING -

250 mL (8 oz.) / 12 per case

### - CATALOG NUMBERS -

<u>Item</u>	<u>Fisher Cat. No.</u>	<u>FoxWest Sales Cat. No.</u>
8 oz.	04-355-135	1409

---

**For Additional Technical Information call  
FoxWest Sales Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)**

---

FoxWest Sales, Inc.

Revised 07/08/09