

**- BRIEF DESCRIPTION -**

*RNase Displace removes RNase, and unwanted DNA. It works instantly so there is no waiting. RNase Displace eliminates the time-consuming method of autoclaving. RNase Displace can be used on all surfaces including glass, plastic, metal, and ceramic. RNase Displace is phosphate-free, safely disposable, non-toxic, and totally rinsable for the decontaminated surface.*

**- USE PROCEDURE -**

**Use RNase Displace full strength.**

**Method 1:** Thoroughly wet item to be decontaminated with RNase Displace - ensure all surfaces are coated. Swirl for a few seconds, drain, and immediately rinse twice with the RNase-free water.

**Method 2:** Soak item in RNase Displace until needed. Drain and rinse as above.

**Method 3:** Soak an RNase-free wipe in RNase Displace, wipe over surface to be decontaminated. Rinse as above.

**For total radioactive decontamination,** soak item in RNase Displace for 2-12 hours. Rinse as above. Wipe over benchtops and sinks to remove residual contamination.

**- CAUTION -**

**Alkaline!** Severe eye irritant, mild skin irritant. Wear eye and skin protection when handling RNase Displace. In case of contact with eyes, immediately flush with water for fifteen minutes and contact a physician immediately.

**- PACKAGING -**

250 mL (8 oz.) / 12 per case

950 mL (32 oz.) / 12 per case

One Gallon / 4 per case

**- CATALOG NUMBERS -**

<u>Item</u>	<u>Fisher Cat. No.</u>	<u>FoxWest Sales Cat. No.</u>
8 oz.	04-355-137	1408
32 oz.	04-355-138	1432
1 gallon	04-355-136	1401

**For Additional Technical Information call  
FoxWest Sales Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)**

FoxWest Sales, Inc.

Revised 07/08/09