



NoCount[®]

**Radioactive Decontaminant for Hands
Technical Data Sheet**

- BRIEF DESCRIPTION -

NoCount aerosol removes radioactive contaminants from hands and skin. NoCount's gentle foam formula helps to lift and trap particles so they can be easily rinsed away with water. Effective for situations where there is accidental contamination of skin.

- USE PROTOCOL -

Shake can well. Dispense No Count onto hands and rub into skin thoroughly making sure foam contacts all contaminated skin surfaces. Rinse with water. Repeat if necessary.

- STORAGE -

Contents under pressure. Do not store near heat source. Do not store above 120° F.

- PRECAUTIONARY STATEMENT -

Eye irritant. Avoid contact with eyes. Always wear gloves when handling radioactive materials.

- PACKAGING -

18 oz. spray can / 12 per case

- CATALOG NUMBER -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	18 oz.	6618

**For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or
See our website at www.deconlabs.com**

[®]Decon Labs, Inc.

Rev. 11/2006