



## Radioactive Decontaminant for Hard Surfaces Technical Data Sheet

### - BRIEF DESCRIPTION -

*NoCount aerosol spray removes radioactive contaminants from benchtops and other hard surfaces. NoCount's foaming action helps to lift and trap particles so they can be easily wiped or rinsed away. Safe to use on all hard surfaces.*

### - USE PROTOCOL -

Shake can well. Spray contaminated area liberally while holding can 12 inches from the contaminated surface. Wait 30 seconds, then wipe away with disposable absorbent material or rinse away with warm water. Repeat if necessary.

### - APPLICATIONS -

For removal of radioactive contaminants from sinks, bench-tops, acrylic shields, glassware, plasticware, electrophoresis gel boxes, sequencing units, etc.

### - STORAGE -

Contents under pressure. Do not store near heat source. Do not store above 120° F.

### - PRECAUTIONARY STATEMENT -

Eye and skin irritant. Wear gloves while handling radioactive materials. Dispose of radioactive waste according to local, state and federal regulations.

### - PACKAGING -

18 oz. spray can / 12 per case

### - CATALOG NUMBER -

<u>Manufacturer</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	18 oz.	6518

**For Additional Technical Information call  
Decon Customer Service  
(800) 332-6647, or  
See our website at [www.deconlabs.com](http://www.deconlabs.com)**

®Decon Labs, Inc.

Revised 11/2006