



Iradecon Wipes

Technical Data Sheet
EPA Reg. No. 69687-1-56753

- BRIEF DESCRIPTION -

Iradecon Wipes are pre-saturated with 0.525 % Sodium hypochlorite solution. Iradecon Wipes can be used to disinfect medical, dental, and laboratory counters as well as any hard, non-porous surface. Each individual, sealed packet contains a pre-moistened towelette, 9 x 9 inches. Iradecon Wipes are registered with the US EPA to meet bacteriocidal, virucidal, and fungicidal efficacy standards.

- USE PROTOCOL -

Open the Iradecon Wipes packet. Remove the pre-moistened, 9x9-inch wipe and unfold. Simply wipe the desired surface to be decontaminated assuring the surface is completely wet. Allow surface to air dry. Discard used Iradecon wipe in a designated waste container. It is a violation of Federal law to use this product in a manner inconsistent with its labeling.

- ACTIVE INGREDIENTS -

0.525 % Sodium hypochlorite solution

- STORAGE AND DISPOSAL -

Store in cool area. Do not reuse wipes. Dispose of used wipe and empty dispenser in trash.

- PRECAUTIONARY STATEMENTS -

**KEEP OUT OF REACH OF
CHILDREN
CAUTION**

FIRST AID: If contact with the eyes occurs, hold eye open and rinse slowly and gently with water for 15-20 minutes. Remove contact lenses. Get prompt medical attention. If swallowed, drink large amounts of water. Do not induce vomiting. Get prompt medical attention. If contact with skin occurs, wash with plenty of soap and water.

- PACKAGING -

100 wipes per package / 8 packages per case

- CATALOG NUMBER -

<u>Manufactured for:</u>	<u>Item</u>	<u>Cat. No.</u>
Decon Labs, Inc.	100 wipes	4600

For Additional Technical Information call
Decon Customer Service
(800) 332-6647, or

See our website at www.deconlabs.com

®Decon Labs, Inc.

Revised 04/01/2004